# **OPERATING INSTRUCTIONS**

for

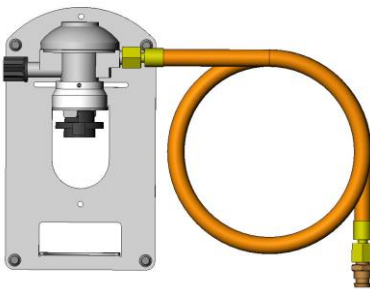
## **Gas Cartridge Adapter CP250**

Part no. 144051


and

## **Gas Cartridge Adapter CV300/470**

Part no. 144052

**CP 250**

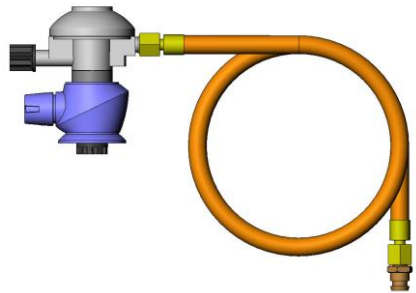**CV 300/470**

## Safety Precautions

### Possible dangers if the safety precautions are not observed

The **Gas Cartridge Adapters CP250 and CV300/470** are state-of-the-art with respect to current technology and safety. Some danger may arise if untrained personnel use them. All persons working with the FIREBOY and the **Gas Cartridge Adapters** must have read and understood the operating and safety instructions or must have been instructed by the supervisors so that safe operation is guaranteed.

### Changes in design or configuration

Changes to the **Gas Cartridge Adapters** are not allowed, neither in design nor in any of their safety aspects. If alterations are made, INTEGRA Biosciences AG will be relieved from all liability for any resulting damage.

## Proper Use

- The described **Gas Cartridge Adapters** are only to be used with the INTEGRA Biosciences AG FIREBOY eco and FIREBOY plus. They are DVGW certified in combination with the FIREBOY only.
- Use only original parts from INTEGRA Biosciences AG for the **Gas Cartridge Adapters** and the specified gas cartridges.
- The **Gas Cartridge Adapters** are only to be used in an environment with temperatures from 10°C to 40°C.
- The **Gas Cartridge Adapters** may not be used without observation. If the **Gas Cartridge Adapters** are not under observation, all valves of the **Gas Cartridge Adapters** must be closed.
- If the **Gas Cartridge Adapters** are not in use, keep them free of dust or dirt (obstruction of connectors).
- Full and empty gas cartridges should be stored in a well-ventilated area.

## Installation of the Gas Cartridge Adapters

1. Prior to connecting the **Gas Cartridge Adapter**, consult the operating instructions of the FIREBOY eco/plus.
2. Check the **Gas Cartridge Adapter** to make sure that the parts don't have visual damages.
3. Turn off FIREBOY.

# INTEGRA

4. Connect the brass connector on the tubing to the gas inlet of the FIREBOY making sure it audibly clicks into position. Check correct positioning by pulling on the tubing.

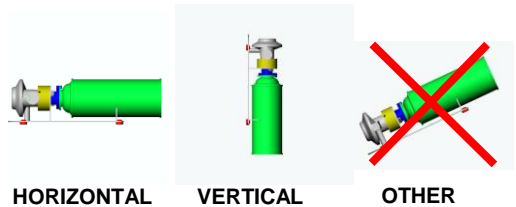
5. Close all valves at the pressure regulator of the **Gas Cartridge Adapter**.

## Mounting the gas cartridge to the Gas Cartridge Adapter CP250

(Part No. 144051)

1. Read the user instruction on the gas cartridge before you connect the gas cartridge to the adapter.
2. Place **Gas Cartridge Adapter** with the gas cartridge in **HORIZONTAL** (work bench) or **VERTICAL** (wall mount) position. Do not move it freely around. The **Gas Cartridge Adapter** must be kept at a distance of minimum 0.5 m from the FIREBOY or any other heat source.

3. Open valve at the pressure regulator of the **Gas Cartridge Adapter**.




Don't open the valves before the tubing is connected to the FIREBOY according to instructions.

4. Turn on the FIREBOY as described in the FIREBOY operating instructions.

## Mounting the gas cartridge to the Gas Cartridge Adapter CV300/470

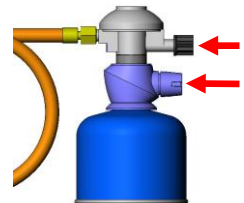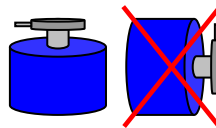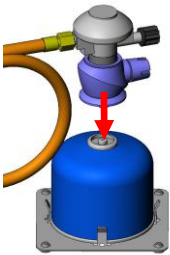(Part No. 144052)

1. Read the user instruction on the gas cartridge before you connect the gas cartridge to the adapter.
2. Use the appropriate gas cartridge stand (CV300 small, CV470 large).
3. Press the **Gas Cartridge Adapter** into the gas cartridge.
4. Gently screw the **Gas Cartridge Adapter** clockwise until you hear a click about one sixth of a turn. Do not tighten the adapter further, as you may damage it.
5. Use the **Gas Cartridge Adapter** only in upright position.
6. Open both valves of the **Gas Cartridge Adapter** by turning counter clockwise as far as the handle.

Don't open the valves before the tubing is connected to the FIREBOY according to instructions.

7. Turn on the FIREBOY according to the FIREBOY operating instructions.

## Removing the gas cartridge from the Gas Cartridge Adapters

1. Turn off the FIREBOY.
2. Close both valves of the **Gas Cartridge Adapter** by turning them clockwise.
3. Hold the gas cartridge and turn the **Gas Cartridge Adapter** counter clockwise about one sixth of a turn. Remove the gas cartridge.

## Removing the Gas Cartridge Adapters from the FIREBOY

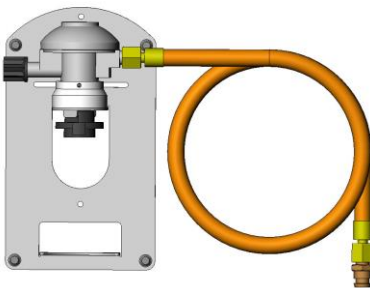
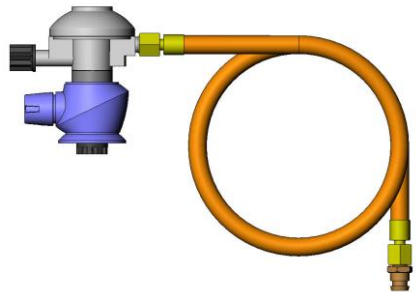
Before removing the **Gas Cartridge Adapter** make sure that all valves of the **Gas Cartridge Adapter** are closed and the gas cartridge is removed.

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