
Bedienungsanleitung ◆ Operating Instructions
Manuel d'utilisation ◆ Manual de instrucciones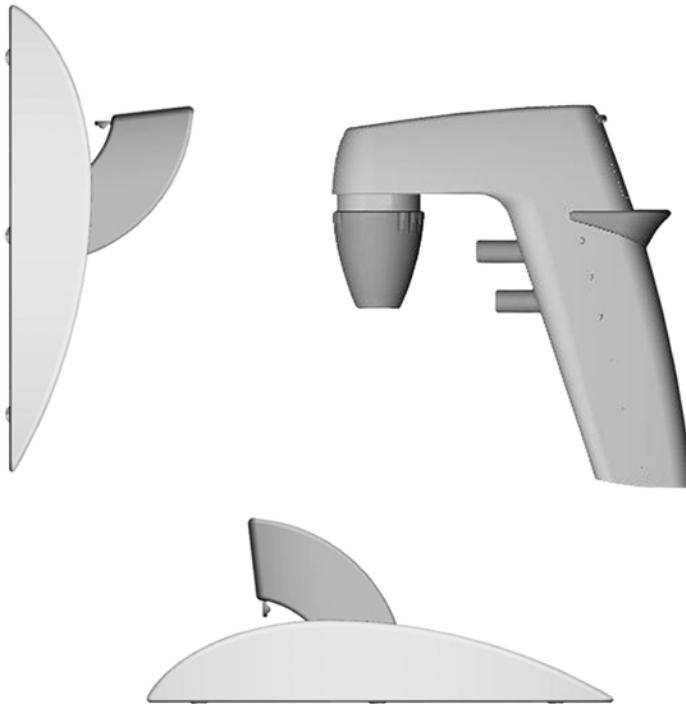

PIPETBOY
comfort

IBS INTEGRA
BIOSCIENCES

Declaration of Conformity

INTEGRA Biosciences (Switzerland) hereby confirms that the product

PIPETBOY comfort

is in accordance with the following directives and standards:

2006/95/EC	Low voltage equipment
2004/108/EC	Electromagnetic compatibility
2002/95/EC	Restriction of Hazardous Substances
2002/96/EC	Waste Electrical and Electronic Equipment
EN 61010-1:2001-02	
EN 61326-1:2006	

CAN/CSA-C22.2 No.1010.1-92

UL Std. No. 3101-1

FCC, Part 15, class A

February 1, 2009

Elmar Morscher
CEO

Thomas Neher
Quality Manager

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to both Part 15 of the FCC Rules and the radio interference regulations of the Canadian Department of Communications. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interface at his own expense.

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Contents

	Page
Safety symbols used	20
Proper use	20
Safety precautions	21
Setting up the PIPETBOY charging station	23
Charging the PIPETBOY	24
Using the PIPETBOY	25
Cleaning/disinfection of the PIPETBOY	26
Replacing the batteries	27
Equipment disposal	28
Troubleshooting	29
Accessories	30
Technical data	31
Manufacturer and service contacts	32

Copyright:

INTEGRA Biosciences

All rights reserved. This manual may not be copied or reproduced in whole or in part in any way without the prior express written permission of INTEGRA Biosciences.

The company reserves the right to make alterations without prior notice.

Notwithstanding the above statements, INTEGRA Biosciences accept no liability for possible mistakes or their consequences arising from this manual.

Safety symbols used

The following are the safety notes and their symbols used in this manual.

Danger

Indicates possible danger of injury or death. Follow the instructions.

Caution

Indicates possible damage to property.

Indication

Gives important information about the correct use of PIPETBOY.

Proper use

The PIPETBOY is designed for the reproducible pipetting and dispensing of liquids. The instrument may not be used for any application other than those specified.

Safety precautions

Possible dangers if the safety precautions are not observed

PIPETBOY is a state-of-the-art instrument with respect to current technology and is safe to use. Some danger may arise if untrained personnel use the instrument. All persons working with the PIPETBOY must read the operating and safety instructions before operating the instrument.

Safety at work

Independently of any safety precautions listed here, all other relevant safety aspects - e.g. GLP, GMP, professional trade associations, Department of Health, trade supervisory authorities - must be observed.

Safety aspects for the user company

The supervisors of operating personnel must have read and understood all the safety precautions involved and must ensure that these are adhered to. They must also ensure that operating personnel also understand and observe all safety rules.

Danger

The instrument must not be used or recharged in an atmosphere with danger of explosion. Highly flammable liquids must not be pipetted (flash point below 0 °C, e.g. Ether, Acetone).

**Caution**

Please dispose the NiMH-batteries according to the relevant provisions.

Restriction of use

The instrument must not be used with solvents whose vapors are incompatible with the materials PA, POM, FPM, NBR, CR, Silicone.

Changes in configuration/design by user

No changes may be made to the PIPETBOY. Any parts that become defective must be replaced by original INTEGRA Biosciences parts.

PIPETBOY may not be altered in design or in any of their safety aspects without the express written permission of INTEGRA Biosciences. In particular, no alterations may be made to any of the protective functions. Any alteration made releases INTEGRA Biosciences from any liability for damage caused.

**Danger**


PIPETBOY and the transformer may not be opened! Repairs must be carried out by INTEGRA Biosciences or a service company authorised by INTEGRA Biosciences.

Setting up the PIPETBOY charging station

Wall mounted


The charging station may be fixed to a wall by the following means:

- double sided self-adhesive tape (1), or,
- magnetic tape attached to the wall (1), or,
- screw through centre of the charging station (2).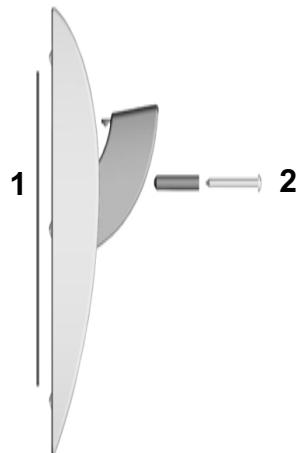

Workbench set-up

Rubber feet on the bottom of the charging station prevent slipping (3).

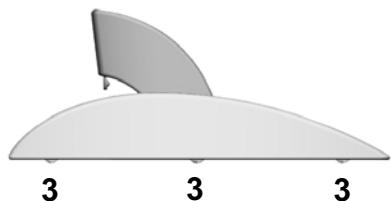Charging the PIPETBOY

- Plug the cable of the mains adapter into the PIPETBOY or the charging station.
- The flashing of the yellow light indicates that batteries are being charged.
- Charge the battery for at least 8 hours before operating the PIPETBOY for the first time. Overcharging is not possible.
- If the yellow light is permanently on, the batteries have discharged to below 10% - recharging is required.
- The PIPETBOY may be used while recharging.


Caution

- Ensure that you are using an original INTEGRA Biosciences mains adapter (100-240 VAC, 50/60 Hz):
US version 156 630
EU version 156 631
UK version 156 632
AU version 156 633
JP version 156 634
- The mains adapter must not be used for any other electrical equipment but the PIPETBOY comfort.

Direct charging of the PIPETBOY

Charging the PIPETBOY in the charging station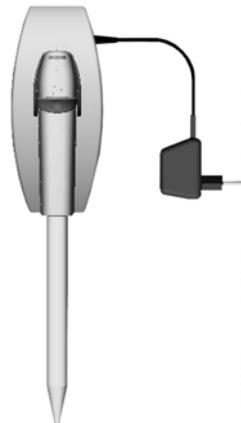

Using the PIPETBOY

Plug in the pipette

- Pipette \geq 2 ml : place the pipette mount (located inside the nose piece) with the small hole on top (1).
- Pipette < 2 ml : place the pipette mount with the large hole on top (2).

To fill the pipette:


- Press the filling control button (3).

To discharge the pipette:

- Press the discharge control button (4).

Regulation of the suction and discharge speed

- By using the pump speed control (5).
- Fine control over the fluid meniscus in the pipette can be achieved by applying varying finger pressure to the operating buttons, as required.

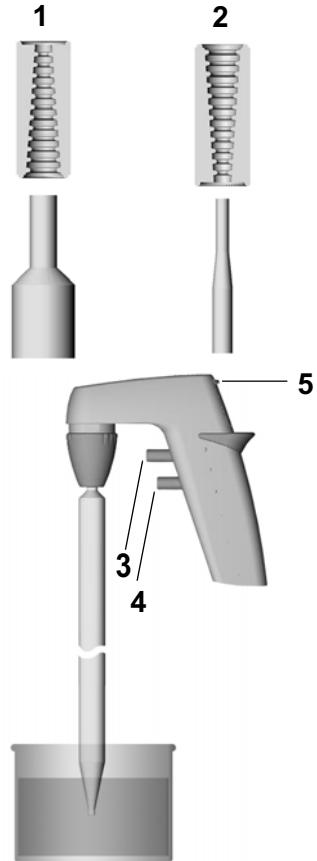Cleaning/disinfection of the PIPETBOY

Danger

Always disconnect the PIPETBOY from the power supply before cleaning.

Caution

Never use acetone to clean the PIPETBOY.


Use only commercially available cleaning agents, such as alcohol (isopropyl alcohol) to clean your PIPETBOY.

Disinfecting the PIPETBOY

Pipette nose piece (1), pipette mount (2) and filter rubber (4) are autoclavable at 121°C.

Changing the filter

The sterile filter (3) should be changed every three months, or if the suction power of PIPETBOY becomes too low.

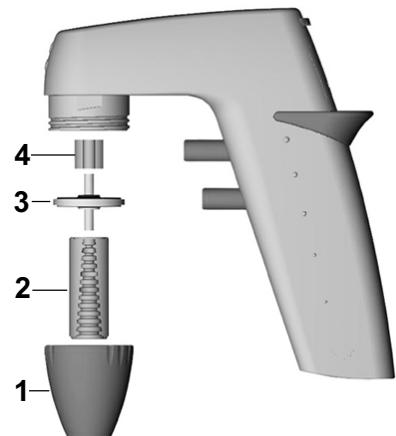Replacing the batteries


Caution

Only use original INTEGRA Biosciences rechargeable batteries (1.2 V AAA NiMH/ 700 mAh).

English

Replacement procedure

- Take out the silicone plug (1).
- Open the screw (2).
- Pull out the flexor (3).
- Open the battery cover (4).
- Change the batteries (1.2 V AAA NiMH / 700 mAh).
- **Check the correct polarity of batteries (+/-) before closing the cover.**

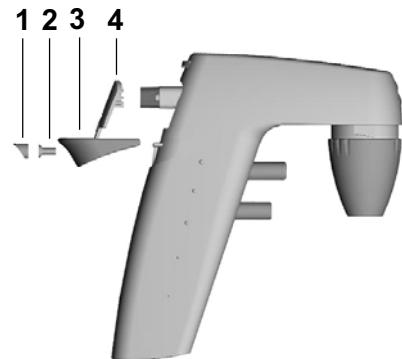Equipment disposal

The packaging of the PIPETBOY *comfort* is labelled with the “crossed-out wheeled bin” symbol to indicate that this equipment must not be disposed of with unsorted municipal waste. Instead, it is your responsibility to correctly dispose of your waste equipment by handing it over to an authorised facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical, and/or radiological contamination so as to protect from health hazards the persons involved in the disposal and recycling of equipment.

For more information about where you can drop off your waste equipment for recycling, please contact your local dealer from whom you originally purchased the product or your local council.

By doing so, you will help conserve natural resources and you will ensure that your waste equipment is recycled in a manner that protects human health and the environment.

Thank you!


Troubleshooting

Problem	Possible cause	Solution
Yellow lamp lights up permanently.	<ul style="list-style-type: none">Batteries are discharged.	<ul style="list-style-type: none">Charge the batteries.
Liquid is not aspirated.	<ul style="list-style-type: none">Missing or defective batteries.	<ul style="list-style-type: none">Check that batteries have been inserted.Charge the batteries.Replace the batteries.
Liquid is not aspirated.	<ul style="list-style-type: none">Pipettor nose piece not tight.	<ul style="list-style-type: none">Tighten the nose piece.
Liquid is not aspirated.	<ul style="list-style-type: none">The filter in the pipettor nose piece is wet and/or clogged.	<ul style="list-style-type: none">Replace the filter.

Accessories

Ordering information

	Art. No.
1. Filter rubber	153 225
2. Filter sterile (0.45 µm)	153 015
Filter not sterile (0.45 µm)	153 016
3. Filter sterile (0.2 µm)	156 608
Filter not sterile (0.2 µm)	156 607
3. Pipette mount, silicone	151 020
4. Pipettor nose piece (classic)	156 100
Pipettor nose piece (volcano)	156 101
Pipettor nose piece (pacific)	156 102
Pipettor nose piece (sahara)	156 103
Pipettor nose piece (royal)	156 104
5. Batteries (2 pieces 1.2 V AAA NiMH/700 mAh)	156 151
6. Battery cover set (classic)	156 110
Battery cover set (volcano)	156 111
Battery cover set (pacific)	156 112
Battery cover set (sahara)	156 113
Battery cover set (royal)	156 114
7. Charging station (classic)	156 050
Charging station (volcano)	156 051
Charging station (pacific)	156 052
Charging station (sahara)	156 053
Charging station (royal)	156 054
8. Mounting kit (magnetic and self-adhesive tape, screw)	156 150
9. Mains adapter (100-240 VAC, 50/60 Hz):	156 630
US version	156 631
EU version	156 632
UK version	156 633
AU version	156 634
JP version	156 634

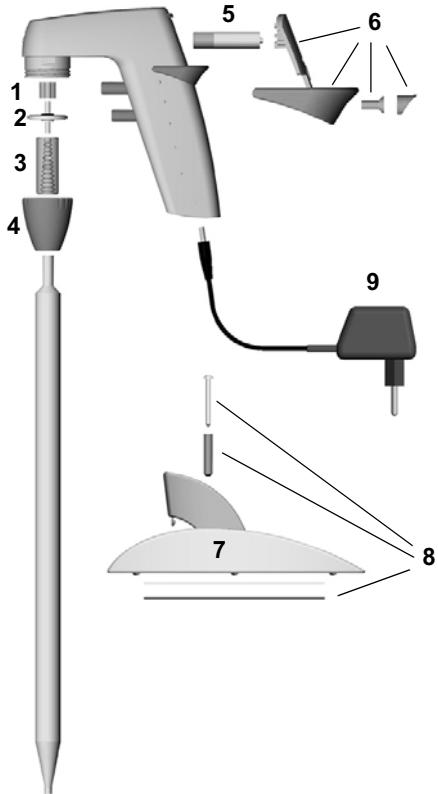Technical data

Dimensions (L x H x W)

PIPETBOY comfort:	130 x 145 x 35 mm
Charging station:	175 x 75 x 80 mm
Mains adapter:	75 x 45 x 45 mm

Weight

PIPETBOY comfort:	190 g
Charging station:	240 g
Mains adapter:	190-300 g

Power supply

Mains adapter:	100-240 VAC, 50/60 Hz (input) 9 VDC, 500 mA (output)
Instrument:	9 VDC or 9 VAC

Battery

Rechargeable batteries 2 x 1.2 V AAA NiMH / 700 mAh

Operating condition

Operation: 5–40°C, max. 80% RH

Storage: -10–50°C, max. 95% RH

UV-Resistance:

If the Pipetbpy will be exposed to ultraviolet light, the plastic housing might discolour and become yellowish. This will not affect the functioning efficiency of the device.

Manufacturer and service contacts

Manufacturer:

INTEGRA Biosciences AG
CH-7205 Zizers
Switzerland
www.integra-biosciences.com
info@integra-biosciences.com

Customer service:

Please contact your local INTEGRA Biosciences AG distributor. You will find their name and address on our web site www.integra-biosciences.com

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