**Bedienungsanleitung • Operator's Manual  
Mode d'emploi • Manuale per l'utilizzatore**

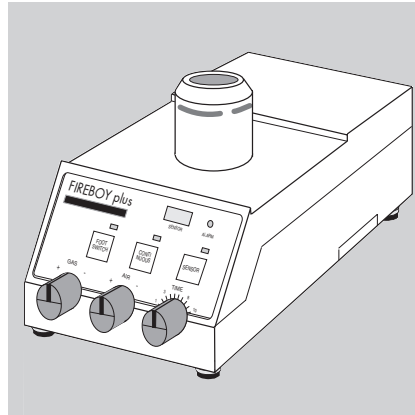
## ***FIREBOY plus***

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**Mobile Bunsen burner with sensor-  
activated ignition**

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**IBS INTEGRA  
BIOSCIENCES**

*Innovative Systems  
for Biotechnology*

**Declaration of Conformity**

**INTEGRA Biosciences** (Switzerland) hereby confirms that the product

**FIREBOY plus**                    **143000**

and

**FIREBOY eco**                    **143100**

is in accordance with the following directives and standards:

**73/23/EEC**      Low voltage equipment

**89/336/EEC**    Electromagnetic compatibility

**2002/95/EC**    Restriction of Hazardous Substances

**2002/96/EC**    Waste Electrical and Electronic Equipment

**IEC/EN 61010-1:2001-02**

**IEC/EN 61326-1+A1:1998+A2:2000**

**CAN/CSA-C22.2 No. 1010.1-92**

**UL Std. No. 3101-1**

**FCC, Part 15, class A**

August 31, 2006

Elmar Morscher  
CEO

Thomas Neher  
Quality Manager

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to both Part 15 of the FCC Rules and the radio interference regulations of the Canadian Department of Communications.

These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications.

Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interface at his own expense.

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## Contens

<b>Delivery schedule .....</b>	<b>2</b>
<b>Appropriate use .....</b>	<b>4</b>
<b>Safety precautions .....</b>	<b>6</b>
<b>Safety features .....</b>	<b>10</b>
<b>Useable types of gas .....</b>	<b>12</b>
<b>Battery .....</b>	<b>14</b>
Charging the battery via a separate charger .....	14
Charging the battery in FIREBOY plus .....	18
<b>Connecting FIREBOY plus to the gas supply .....</b>	<b>18</b>
Selecting the burner nozzle .....	18
Connecting to the central gas supply or to a gas cylinder .....	20
Gas feed via gas cartridge .....	22
Changing the gas cartridge .....	24
Switching from cartridge operation to permanent gas connection .....	26
Setting the right flammable mixture .....	28
<b>FIREBOY plus – operating modes .....</b>	<b>28</b>
Operating mode "FOOT SWITCH" .....	28
Operating mode "CONTINUOUS" .....	30
Operating mode "SENSOR" .....	32
<b>Working with liquids .....</b>	<b>36</b>
<b>Equipment disposal .....</b>	<b>36</b>
<b>Troubleshooting .....</b>	<b>40</b>
<b>Operation instructions FIREBOY eco .....</b>	<b>42</b>
Connecting to the mains .....	42
Connecting the foot switch .....	44
Preparing the instrument for operation .....	44
<b>Technical Data .....</b>	<b>46</b>
<b>Accessories .....</b>	<b>48</b>

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## Delivery schedule

Check all components against delivery schedule and ensure that they are complete and undamaged.

### **FIREBOY plus:**

- 1 FIREBOY plus instrument
- 1 Power supply
- 1 Nozzle P60 propane/butane gas (fitted ex-factory)
- 1 Nozzle N80 natural gas
- 1 Nozzle M190 mixed gas
- 1 Battery rechargeable
- 1 Tube adapter Ø 10 mm
- 1 Hose clip
- 1 User manual

### **FIREBOY eco:**

- 1 FIREBOY eco instrument
- 1 Power supply
- 1 Nozzle P60 propane/butane gas (fitted ex-factory)
- 1 Nozzle N80 natural gas
- 1 Nozzle M190 mixed gas
- 1 Foot switch
- 1 Tube adapter, Ø 10 mm
- 1 Hose clip
- 1 User manual

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## *Appropriate use*

FIREBOY plus/eco is an automatic Bunsen burner for flame-sterilising laboratory materials made of glass or metal. Never use for any other purpose!

### **Changes in Design or Configuration**

No changes may be made to FIREBOY. Original INTEGRA Biosciences parts must be used to replace any parts that become defective.

FIREBOY may not be altered in design or in any of its safety aspects. If alterations are made, INTEGRA Biosciences will be relieved from all liability for any resulting damage.

**FIREBOY must not be opened! Repairs must be carried out by INTEGRA Biosciences or by a service company authorized by INTEGRA Biosciences.**

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**ATTENTION: Hot**

**Caution in the "sensor" mode**

- The equipment may be operated only under constant supervision.
- Always wear protective spectacles when working with FIREBOY plus/eco.
- Use only approved and tested gas tubing.
- Use only the nozzles supplied with the instrument.
- Do not use near flammable material or in highly combustible areas.
- Never remove the gas cartridge adapter whilst it is connected to the gas cartridge.

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- Only use original INTEGRA Biosciences AC adapter.
- Only use original INTEGRA Biosciences rechargeables batteries.  
Never use non-rechargeables batteries!
- Please dispose the battery and the gas cartridge according to the relevant provisions.

**The gas supply must be closed off when:**

- the instrument is being transported or installed
- the timer is being installed at the base of the instrument
- the gas cartridge is being changed
- work has been completed!
- the battery is being changed
- the adapter is being changed

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## Safety Features

- The burner chamber and valve housing are monitored by a temperature sensor.
- If the temperature exceeds the safety level, the gas valve is automatically closed off.
- FIREBOY plus also has a flame monitor: if the flame has not ignited 3 seconds after switching on, the gas valve is automatically closed off.
- If the flame is inadvertently extinguished, the gas valve is closed off automatically within 3 seconds.

### Safety devices for "Sensor" and "Foot switch" operating modes

- Both functions switch off automatically 5 minutes after last use.

### Safety device for operating mode "Continuous"

- The flame is switched off automatically after 10-60 minutes, depending on the timer setting.

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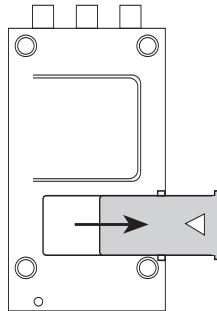
## Battery

FIREBOY plus can be operated independently of the power supply by means of the integrated battery. The battery can be recharged either separately with a battery charger or in FIREBOY plus.

**The battery should be charged for 14 hours on being charged for the first time. All subsequent charging takes place via a rapid charger and is complete in approximately 3 hours.**

### Charging the battery via a separate charger

- Slide the cover of the battery compartment in the direction of the arrow.

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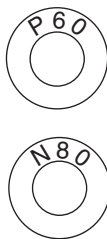
## Charging the battery in FIREBOY plus

FIREBOY plus incorporates a charge protection device to prevent the battery from being overcharged. This guarantees maximum battery life.

FIREBOY plus can also be operated without a battery via the power supply.

- Connect the charger plug to socket ① on FIREBOY plus and then connect to the mains.

## *Connecting FIREBOY plus to the gas supply*

### Selecting the burner nozzle

Accordingly the correct nozzle must be used. See chapter "Useable types of gas".

The nozzle for propane/butane gas is fitted ex-factory.

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
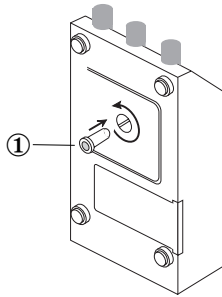
- Remove screw ① from the base of the instrument.
- Remove the nozzle.
- Insert the required nozzle.
- Replace screw ① tightly in position.

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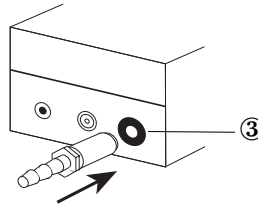
## Connecting to the central gas supply or to a gas cylinder

- Connect the gas adapter to gas inlet ③, making sure it audibly clicks into position.
- Connect the pipe from the central supply or from the gas cylinder to the adapter and fix with a clip if necessary.

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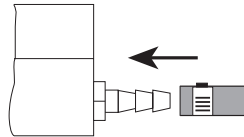### Changing the adapter:

**Switch off all operating modes - all green LEDs must be extinguished !**

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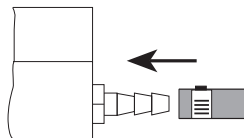
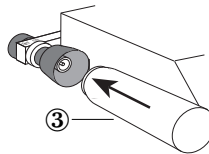
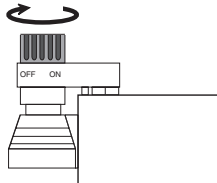
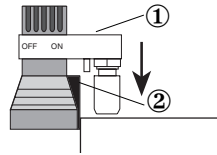
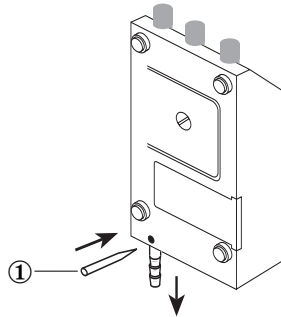
- Use a pointed tool ① to press the adapter mechanism inwards and remove the adapter by pulling towards the rear.
- Insert the requested new adapter until it audibly clicks into position.

## Gas feed via gas cartridge

When using a gas cartridge, make sure that nozzle P 60 has been fitted.

- Insert the gas cartridge adapter ① into the Gas Inlet at the rear of the instrument until it audibly clicks into position. Ensure that the smooth side of cap ② is flush with the FIREBOY plus housing.
- Turn the milled knob on the cartridge adapter fully clockwise in the "OFF" direction.
- Insert the gas cartridge ③ into the black plastic part of the adapter until it also audibly clicks into position.

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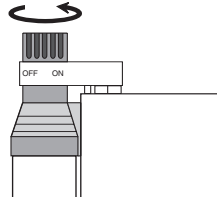
### Then open the gas cartridge adapter:

- Turn the milled knob on the cartridge adapter fully anticlockwise in the "ON" direction.

The gas cartridge is now open and ready for use.

The flame can now be ignited using any of the three operating modes.

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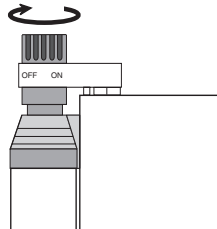
## Changing the gas cartridge

### Caution:

#### Before proceeding, close off the gas cartridge adapter:

- Turn the milled knob on the cartridge adapter fully clockwise in the "OFF" direction.
- Remove the empty cartridge.
- Fit a new cartridge.
- Turn the milled knob on the cartridge adapter fully anticlockwise in the "ON" direction.

The gas cartridge is now ready for use (**Caution: gas will begin to flow**).

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## Switching from cartridge operation to permanent gas connection

### Before proceeding, close off the gas cartridge adapter:

- Turn the milled knob on the cartridge adapter fully clockwise in the "OFF" direction.
- Remove the gas cartridge.

### Then remove the gas cartridge adapter:

- Using a pointed tool, press locking mechanism ① on the underside of the FIREBOY plus base and remove the adapter by pulling to the rear.

- Fit the adapter appropriate for the laboratory gas supply to the gas inlet (see page 22). If necessary, change the nozzle.

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## Setting the right flammable mixture

The "GAS" and "AIR" knobs can be rotated to set the right gas/air mixture in conjunction with observation of the flame itself.

The "GAS" knob can also be used to adapt to the particular type of gas being used.

## *FIREBOY plus – operating modes*

FIREBOY plus can be operated in three modes:

- "FOOT SWITCH"
- "CONTINUOUS"
- "SENSOR"

### Operating mode "FOOT SWITCH"

In this operating mode, the flame is ignited only when the foot switch is activated. The flame burns only as long as the foot switch is depressed.

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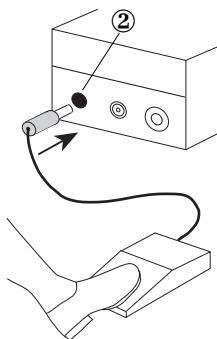
In the "FOOT SWITCH" mode, the flame is extinguished automatically 5 minutes after the last foot switch activation. The green LED is also extinguished.

#### **Preparing for the "FOOT SWITCH" operating mode:**

- Connect the foot switch plug to socket ② at the rear of the instrument.
- Press the "FOOT SWITCH" key on the keyboard. The green LED will light.

The flame can now be ignited and extinguished by activating the foot switch.

#### **Switching off in the "FOOT SWITCH" mode:**

- Press the "FOOT SWITCH" key. The green LED will extinguish.

#### **Operating mode "CONTINUOUS"**

In this operating mode, the duration of flame operation can be regulated by a timer that can be set from 10 to 60 minutes.

The timer is located under the base of the instrument.

The timer is adjusted ex-factory to approximately 10 minutes.

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### **Setting the flame duration by timer in the "CONTINUOUS" mode:**

**NB: First switch off the gas supply! And switch off all operating modes.**

- Using a screwdriver:
  - Turn the timer to the left (for shorter duration)
  - Turn the timer to the right (for longer duration).

### **Switching on in "CONTINUOUS" mode:**

- Press the "CONTINUOUS" key on the keyboard. The green LED will light. FIREBOY will ignite and continue burning until the set period has expired.

### **Switching off in "CONTINUOUS" mode:**

- Press the "CONTINUOUS" key on the keyboard. The flame and the green LED will extinguish.

### **Operating mode "SENSOR"**

In this operating mode, the flame will ignite when an object slowly passes the sensor window (infrared beam) at a maximum distance of 6-8 cm.

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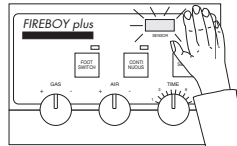
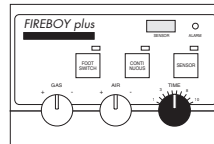
The flame duration can be set between 3 and 13 seconds with the rotary knob (TIME) on the keyboard.

The "SENSOR" operating mode remains active for 5 minutes after the last ignition; if the flame is not re-ignited after this time, the "SENSOR" mode will be switched off and the green LED above sensor will be extinguished.

### Switching on in "SENSOR" mode:

- Press the "SENSOR" key on the keyboard. The green LED will light.
- Set the flame duration between 3 - 13 seconds with the "TIME" knob.

- Using a hand or appropriate object, pass the sensor window. The flame will ignite and continue burning until the set period has expired.

### Switching off in "SENSOR" mode:

- Press "SENSOR" on the keyboard. The green LED will extinguish.

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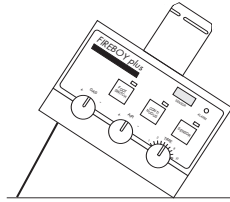
## Working with liquids

The folding stand at the base of the instrument can be used to incline the instrument. This protects the burning chamber from contamination when working with liquids.

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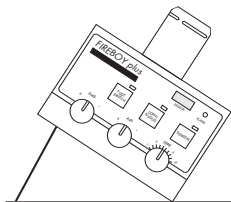
## Equipment disposal

The packaging of the FIREBOY is labelled with the "crossed-out wheeled bin" symbol to indicate that this equipment must not be disposed of with unsorted municipal waste. Instead, it is your responsibility to correctly dispose of your waste equipment by handing it over to an authorised facility for separate collection and recycling.

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It is also your responsibility to decontaminate the equipment in case of biological, chemical, and/or radiological contamination so as to protect from health hazards the persons involved in the disposal and recycling of equipment.

For more information about where you can drop off your waste equipment for recycling, please contact your local dealer from whom you originally purchased the product or your local council.

By doing so, you will help conserve natural resources and you will ensure that your waste equipment is recycled in a manner that protects human health and the environment. Thank you!

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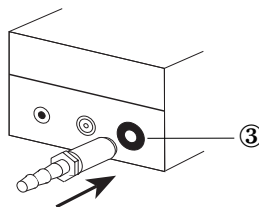
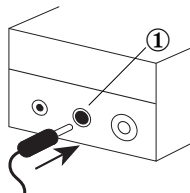
## Operating instructions FIREBOY eco

FIREBOY eco is a mains-dependent instrument and the flame remains ignited as long as the foot switch is activated.

FIREBOY eco has neither a flame monitor nor an overheating monitor.

### Connecting FIREBOY eco to the mains

- Connect the power supply plug to socket ① at the rear of the instrument and connect it to the mains.
- Check that the appropriate nozzle is fitted for the gas in use.
- Connect the gas adapter to gas inlet ③, making sure it audibly clicks into position.

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INTEGRA Biosciences AG