Operation


The vacuum is controlled by the pressure sensitive turquoise control button on the top of the hand piece. By means of the plastic screw opposite the control button, the residual suction level can be set, which will prevent dripping of waste liquids from the tip or pipette. The aspirator is factory pre-set to allow a low level of constant suction; by turning the screw one full turn clockwise, the valve is shut, and residual vacuum is stopped.

Cleaning and servicing

The Hand Operator may be decontaminated with H₂O₂ gas (max. 35 %) for 60 min and is entirely autoclavable. When working with biohazardous material, disinfect the VACUBOY with a 70 % solution of isopropanol or ethanol or a suitable disinfectant before disassembling for autoclaving. Please refer to the explosion drawing.

To clean, immerse the pieces in mild laboratory detergent or 70 % ethanol, but do not use abrasives, as a roughened surface may affect the smooth operation of the control valve. Apply a thin coat of grease MOLYKOTE as needed to the control valve to keep the action smooth (arrow).

When working with salt-rich solutions, rinse out the aspirator hand piece with distilled water after use to avoid build-up of dried deposits.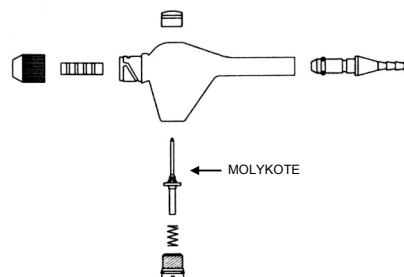

Chemical compatibility

Avoid prolonged contact (days or weeks) with:

- concentrated inorganic acids
- sodium hypochlorite
- methylene chloride
- concentrated aldehydes
- other halogenated hydrocarbons


Resistant to:

- DMSO
- ethanol/methanol
- dimethylformamide
- acetone

Note:

- Flushing the aspirator with ethanol immediately after contact with organic solvents is the best safeguard in all cases.
- For O-ring replacement, use INTEGRA O-rings only

Accessories and consumables

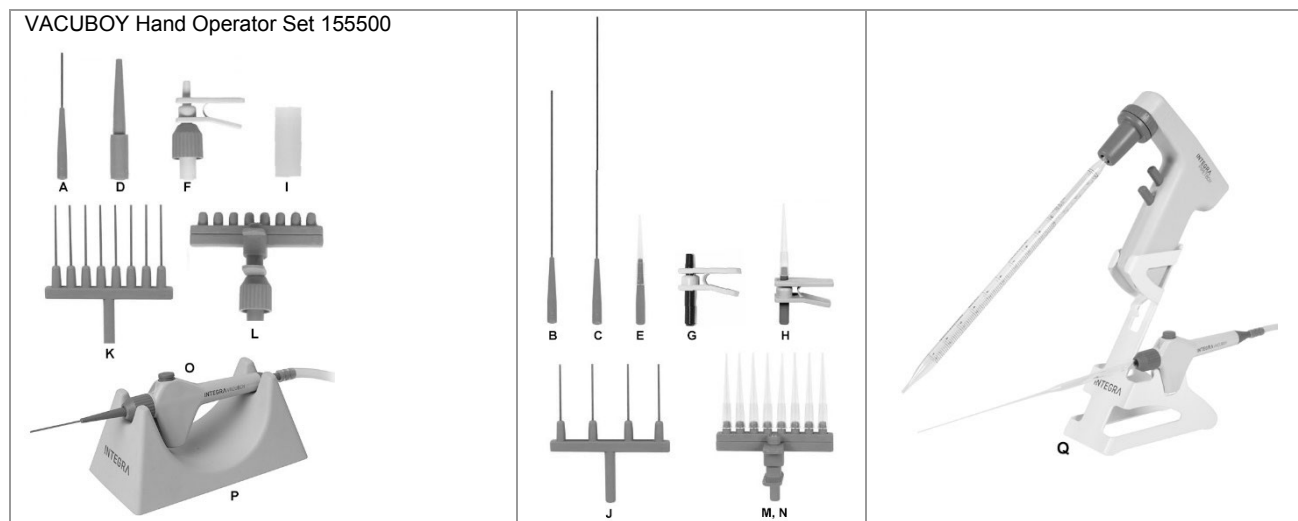VACUBOY Hand Operator: hand operator (O), two aspiration adapters (A and F), rubber adapter for Pasteur pipettes (I), grease (Dow Corning MOLYKOTE HP-870)	155510
VACUBOY Hand Operator Set: hand operator (O), aspiration adapter D, A, K, L, F, rubber adapter for Pasteur pipettes (I), stand for hand operator (P), silicone tubing 1800 mm, grease	155500
A 1-channel stainless steel aspiration adapter 40 mm	155502
B 1-channel stainless steel aspiration adapter 150 mm	155522
C 1-channel stainless steel aspiration adapter 280 mm	155525
D 1-channel aspiration adapter for disposable tips/GripTips (with locking)	155504
E 1-channel aspiration adapter for disposable tips/GripTips (pack of 5)	159023
F 1-channel aspiration adapter with ejector for disposable tips (with locking)	155526
G 1-channel aspiration adapter with ejector for disposable tips	159026
H 1-channel aspiration adapter with ejector GripTips	159027
I Rubber adapter for Pasteur pipettes	155505
J 4-channel stainless steel aspiration adapter 40 mm	155524
K 8-channel stainless steel aspiration adapter 40 mm	155503
L 8-channel aspiration adapter with ejector for disposable tips (with locking)	155520
M 8-channel aspiration adapter with ejector for disposable tips	159024
N 8-channel aspiration adapter with ejector for GripTips	159025
P Stand for VACUBOY hand operator	155501
Silicone tubing 1800 mm	158332
Q Stand for PIPETBOY (acu or pro) and VACUBOY	155065

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