Operation


The vacuum is controlled by the pressure sensitive turquoise control button on the top of the hand piece. By means of the plastic screw opposite the control button, the residual suction level can be set, which will prevent dripping of waste liquids from the tip or pipette. The aspirator is factory pre-set to allow a low level of constant suction; by turning the screw one full turn clockwise, the valve is shut, and residual vacuum is stopped.

Cleaning and servicing

The Hand Operator may be decontaminated with H₂O₂ gas (max. 35 %) for 60 min and is entirely autoclavable. When working with biohazardous material, disinfect the VACUBOY with a 70 % solution of isopropanol or ethanol or a suitable disinfectant before disassembling for autoclaving. Please refer to the explosion drawing.

To clean, immerse the pieces in mild laboratory detergent or 70 % ethanol, but do not use abrasives, as a roughened surface may affect the smooth operation of the control valve. Apply a thin coat of grease MOLYKOTE as needed to the control valve to keep the action smooth (arrow).

When working with salt-rich solutions, rinse out the aspirator hand piece with distilled water after use to avoid build-up of dried deposits.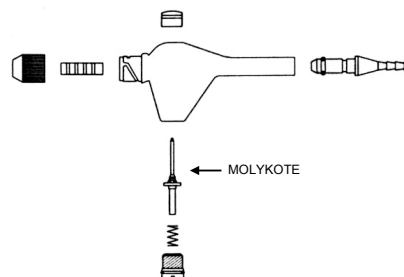

Chemical compatibility

Avoid prolonged contact (days or weeks) with:

- concentrated inorganic acids
- sodium hypochlorite
- methylene chloride
- concentrated aldehydes
- other halogenated hydrocarbons


Resistant to:

- DMSO
- ethanol/methanol
- dimethylformamide
- acetone

Note:

- Flushing the aspirator with ethanol immediately after contact with organic solvents is the best safeguard in all cases.
- For O-ring replacement, use INTEGRA O-rings only

Accessories and consumables

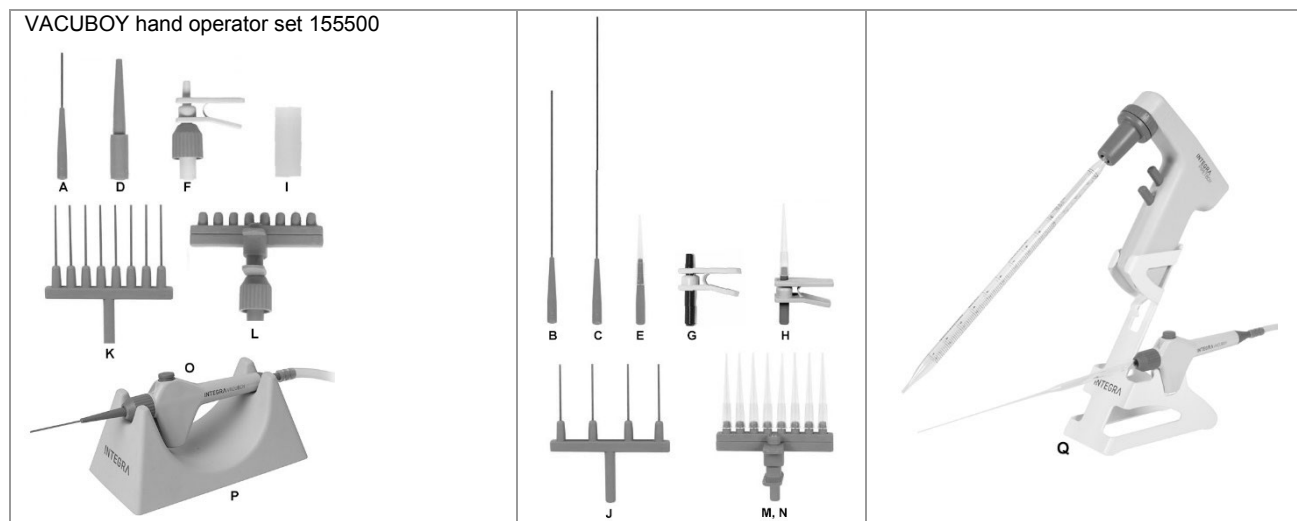| | | |
|--|---|--------|
| VACUBOY hand operator set small: hand operator (O), two aspiration adapters (A and F), rubber adapter for Pasteur pipettes (I), grease (Dow Corning MOLYKOTE HP-870) | | 155510 |
| VACUBOY hand operator set large: hand operator (O), aspiration adapter D, A, K, L, F, rubber adapter for Pasteur pipettes (I), stand for hand operator (P), silicone tubing 1.8 m, grease | | 155500 |
| A | Adapter 1 channel, 40 mm, stainless steel | 155502 |
| B | Adapter 1 channel, 150 mm, stainless steel | 155522 |
| C | Adapter 1 channel, 280 mm, stainless steel | 155525 |
| D | Adapter 1 channel, with locking for tips/GRIPTIPS, plastic | 155504 |
| E | Adapter 1 channel, for tips/GRIPTIPS, plastic, pack of 5 | 159023 |
| F | Adapter 1 channel, with ejector and locking for tips, plastic | 155526 |
| G | Adapter 1 channel, with ejector for tips, plastic | 159026 |
| H | Adapter 1 channel, with ejector for GRIPTIPS, plastic | 159027 |
| I | Adapter for Pasteur pipettes, rubber | 155505 |
| J | Adapter 4 channel, 40 mm, stainless steel | 155524 |
| K | Adapter 8 channel, 40 mm, stainless steel | 155503 |
| L | Adapter 8 channel, with ejector and locking for tips, plastic | 155520 |
| M | Adapter 8 channel, with ejector for tips, plastic | 159024 |
| N | Adapter 8 channel, with ejector for GRIPTIPS, plastic | 159025 |
| P | Stand for VACUBOY hand operator | 155501 |
| | Silicone tubing, Ø 8 mm, 1.8 m, VACUBOY to bottle | 158332 |
| Q | Stand for PIPETBOY (acu or pro) and VACUBOY | 155065 |

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