

Automated qPCR master mix preparation with the D-ONE single channel pipetting module and the ASSIST PLUS pipetting robot

Introduction

Automating qPCR master mix preparation is often considered extremely tedious, due to the need to constantly change program set-up based on varying master mix volumes and sample numbers. However, the D-ONE single channel pipetting module for the ASSIST PLUS pipetting robot simplifies the programming of automated qPCR master mix preparation protocols down to a single step. The D-ONE individually reaches tubes of different qPCR master mix components, while the VIALAB program takes care of all calculations and provides automatic tip selection.

This solution allows master mixes to be pipetted into any tube or reservoir, enabling large-scale qPCR master mix preparation, with [subsequent distribution to plates with multichannel pipettes](#) to speed up the process. Here, we verify a SYBR® Green qPCR protocol prepared directly in a divided reservoir using the D-ONE on the ASSIST PLUS.

Key benefits:

- Fully automated liquid handling of various qPCR master mix components on the ASSIST PLUS pipetting robot with the D-ONE single channel pipetting module ensures the reproducibility of results.
- qPCR master mix components and volumes can be easily adjusted in VIALAB to perform different qPCR protocols. The final calculation is automatically done by VIALAB according to the selected reaction count.
- The D-ONE's liquid level detection allows the use of reagents with different aliquot volumes, and informs the user if there is insufficient liquid. Furthermore, the automated tip selection guarantees high precision when pipetting different volumes.
- qPCR master mixes can be pipetted into various target vessels, including INTEGRA's multichannel reservoirs with a very low dead volume, to speed up the plating of master mixes using a multichannel pipette.

Overview: How to automate protocols for SYBR Green qPCR master mix preparation

The single channel 5-1250 µl D-ONE single channel pipetting module with 125 µl and 1250 µl sterile, filter, GRIPTIPS is mounted on the ASSIST PLUS pipetting robot.



An INTEGRA divided polypropylene reagent reservoir with SureFlo™ anti-sealing array is placed onto the dual reservoir adapter on deck position A. An INTEGRA tube rack with nuclease-free water (A1), SYBR Green master mix reagent (B1), forward primer (C1) and reverse primer (D1) is placed on position B (**Figure 1**).

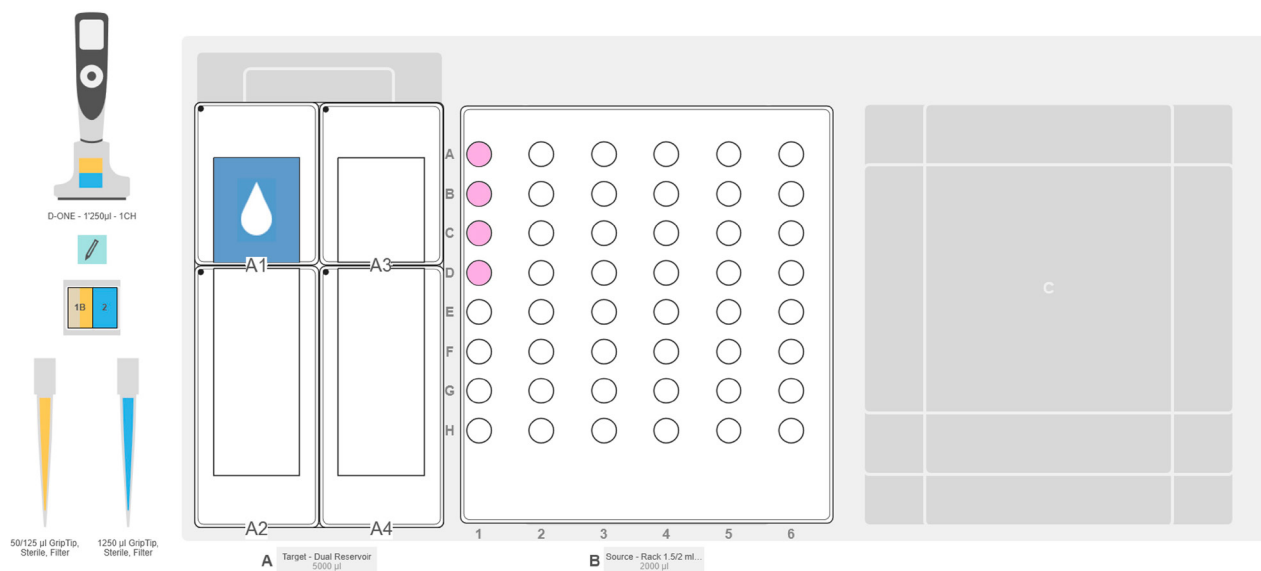


Figure 1: Deck set-up for qPCR master mix preparation in divided reagent reservoir. **Position A:** Target – INTEGRA dual reservoir adapter with divided reagent reservoir (blue). **Position B:** Source – INTEGRA tube rack for 2 ml tubes with screw caps and 1.5 ml microcentrifuge tubes (pink). **Position C:** empty.

Step by step procedure:

1. Calculation of qPCR master mix components

STEP: Adjusting the number of reactions in VIALAB.

The reaction count is predefined to 98 reactions to fill a full 96 well qPCR plate (9 µl reaction volume, 10 µl dead volume). The user defines the reagent ID and its volume per reaction in VIALAB, which can be changed according to the specific qPCR protocol. By adjusting the reaction count, the transfer volume of all qPCR master mix components is calculated automatically, as well as the selection of the appropriate GRIPTIP (**Figure 2**).

TIPS:

- Once a program with a master mix step is set up, it can be reopened and only the number of reactions needs to be changed prior to the start of the run. No reprogramming is needed, saving time for the operator.
- Depending on the final volume calculated by VIALAB, the tip type (125/1250 µl) is automatically chosen, further alleviating the need for user input.

Enter the number of reactions

Reactions		
-	98	+

Reagent ID	Source		Volume (1*) [µl]	Reactions	Validation
	Pos.	Well			
				98	
Nuclease-free Water	B	A1	3	294	✓
SYBR Green Master Mix	B	B1	5	490	✓
Forward Primer	B	C1	0.5	49	✓
Reverse Primer	B	D1	0.5	49	✓
Target	A	A1	9	882 µl	

Figure 2: In the master mix step, VIALAB calculates the volume of qPCR master mix components according to the reaction count.

2. Transfer of qPCR master mix components

STEP: The different components of qPCR master mix are added one by one.

HOW TO: Select and run the VIALAB program 'Master_mix_prep_in_reservoir'. The 5-1250 µl D-ONE pipetting module with 1250 µl sterile, filter GRIPTIPS aspirates 294 µl of nuclease-free water from a 2 ml tube (Position B - A1), and dispenses it into the 5 ml compartment of the divided reservoir Position A (A1). The D-ONE module then changes tips automatically between different reagents, and transfers 490 µl of SYBR Green master mix from B1 to A1 of the divided reservoir. With 12.5 µl sterile, filter GRIPTIPS, the D-ONE then transfers 49 µl each of the forward primer (C1) and reverse primer (D1) into A1 of the divided reservoir. With a new 1250 µl sterile, filter GRIPTIP, the D-ONE mixes 600 µl of SYBR Green master mix three times, in three different offsets within the 5 ml compartment of the divided reservoir (A1).

TIP:

- Different mixing offsets in the reagent reservoir improve the homogeneity of the qPCR master mix components. Mixing volume can be adjusted easily when reducing the reaction count, while mixing cycles can be increased to produce even higher volumes of qPCR master mix.

Results

We used a SYBR Green qPCR protocol to amplify a 250 bp fragment from the fifth variable region of bacterial 16S rRNA gene with the QuantStudio™ 3 (Thermo Fisher Scientific) to demonstrate proper homogeneity when performing qPCR master mix preparation in divided reagent reservoirs. The master mix was prepared both manually in tubes and with the ASSIST PLUS pipetting robot and the D-ONE single channel pipetting module in the divided reservoir. Both qPCR master mix preparations show valid amplification, demonstrating the proper homogeneity of qPCR master mix components.

Figure 3 represents the amplification of a section of six replicates from a qPCR plate using SYBR Green master mix produced with the ASSIST PLUS and the D-ONE in a divided reservoir (green), compared to six replicates using master mix produced manually in tubes (orange). Both qPCR master mix preparations show valid amplification, demonstrating the proper homogeneity of qPCR master mix components.

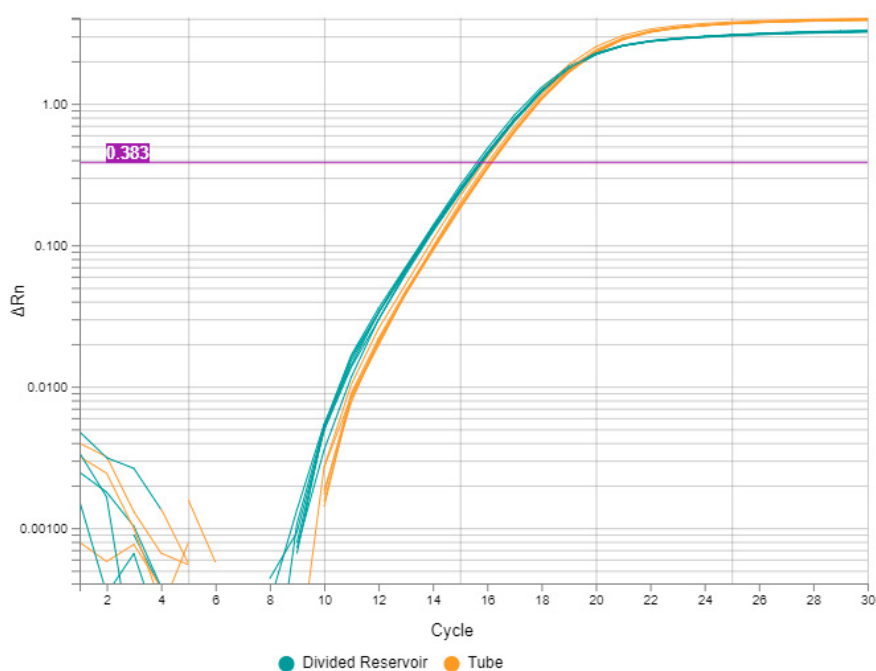


Figure 3: Amplification plot of a 250 bp fragment of the fifth variable region of bacterial 16S rRNA showing a section of six replicates.

Remarks

- **VIALAB software:** VIALAB programs can be easily adapted to your specific labware and protocols, such as when partial plates are needed.
- **Partial plates:** Programs can be adapted at any time to varying sample numbers, giving laboratories total flexibility to meet current and future demands.

Conclusion

- The ASSIST PLUS pipetting robot and D-ONE single channel pipetting module, together with the VIALAB master mix step enables the easy and reliable automation of qPCR master mix preparation, including automatic volume calculations, tip selection and liquid level detection.
- The ASSIST PLUS is the ideal solution for automating qPCR master mix preparation, set-up, and sample addition. It maximizes hands-free time and increases reproducibility by eliminating any variability introduced by operators.
- With its small footprint, the ASSIST PLUS fits into biosafety cabinets to ensure sterile conditions while pipetting precious qPCR master mix components.
- The D-ONE can individually transfer various qPCR master mix components, and effectively mix them in divided reservoirs for large scale qPCR master mix preparation.
- The homogeneity of qPCR master mix components can be guaranteed by adjusting mixing volumes and cycles, setting different mixing offsets, and the unique surface treatment of the SureFlo polypropylene divided reagent reservoir.
- The divided reagent reservoir can then be easily accessed using an 8 channel 12.5 µl VOYAGER adjustable tip spacing pipette to quickly set up qPCR plates.

Materials

Manufacturer	Part Number	Description	Link
INTEGRA Biosciences	4505	ASSIST PLUS base unit	https://www.integra-biosciences.com/global/en/pipetting-robots/d-one-for-assist-plus
INTEGRA Biosciences	4532	D-ONE pipetting module	https://www.integra-biosciences.com/global/en/pipetting-robots/d-one-for-assist-plus
INTEGRA Biosciences	4535	D-ONE Tip Deck	https://www.integra-biosciences.com/global/en/pipetting-robots/d-one-for-assist-plus
INTEGRA Biosciences	4540	Rack for 1.5/2 ml microcentrifuge tubes	https://www.integra-biosciences.com/global/en/pipetting-robots/assist-plus
INTEGRA Biosciences	4547	Dual Reservoir Adapter	https://www.integra-biosciences.com/global/en/pipetting-robots/assist-plus
INTEGRA Biosciences	4356	25 ml Divided Reservoir, Polypropylene	https://www.integra-biosciences.com/global/en/reagent-reservoirs/multichannel-reagent-reservoirs
INTEGRA Biosciences	6465	125 µl Sterile, Filter GRIPTIPS	https://www.integra-biosciences.com/global/en/pipette-tips/griptip-selector-guide
INTEGRA Biosciences	6445	1250 µl Sterile, Filter GRIPTIPS	https://www.integra-biosciences.com/global/en/pipette-tips/griptip-selector-guide

INTEGRA Biosciences AG
7205 Zizers, Switzerland
T +41 81 286 95 30
F +41 81 286 95 33
info@integra-biosciences.com

INTEGRA Biosciences Corp.
Hudson, NH 03051, USA
T +1 603 578 5800
F +1 603 577 5529
info-us@integra-biosciences.com

INTEGRA Biosciences Deutschland GmbH
35444 Biebertal, Deutschland
T +49 6409 81 999 15
F +49 6409 81 999 68
info-de@integra-biosciences.com

INTEGRA Biosciences SAS
95062 Cergy-Pontoise Cedex 1, France
T +33 (0)1 34 30 76 76
F +33 (0)1 34 30 76 79
info-fr@integra-biosciences.com

INTEGRA Biosciences Ltd.
Thatcham, Berks RG19 4EP, UK
T: +44 1635 797000
F: +44 1635 797001
info-uk@integra-biosciences.com

INTEGRA Biosciences Nordic ApS
Vallensbækvej 22A 3TV
Brøndby 2605, Denmark
T + 45 3173 5373
info-nordic@integra-biosciences.com

INTEGRA Biosciences (Shanghai) Co., Ltd.
中国上海自由贸易试验区环科路515号1110室
邮编: 201315
电话: +86 21 5844 7203
info-cn@integra-biosciences.com

インテグラ・バイオサイエンス株式会社
〒101-0031
東京都千代田区東神田 1-5-6
東神田MK第五ビル 3階
T 03-5962-4936
F 03-5962-4937
info-jp@integra-biosciences.com